

Quick Start Guide

MVE Condition Monitoring



Note: Initial registration of logger(s) may take up to 72 hours. Customer Service will contact you once the registration is completed.

Questions? Contact Customer Service at monitor@mvebio.com

CryoBeacon



Temperature



Humidity



Scan the “Register Me” QR code located on the CryoBeacon. If you are unable to scan the code, you may register your logger by visiting <https://register.mvebio.com> and entering the serial number found on the top of the logger under the “Register Me” label.



Download the ‘Tec4App’ available on both the App Store and Google Play Store.



Scan the QR code on the CryoBeacon with Tec4App and read the data immediately.

Once registration is completed, you will receive your login information for the MVECloud.

Quick Start Guide

MVE Condition Monitoring



SmartTag



Temperature



Orientation



Shock




Light



Geo-Position

Display Info

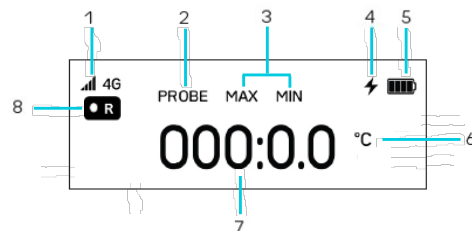

- | | |
|-----------------------|----------------------|
| 1. Connection Icon | 5. Battery Status |
| 2. Probe Indicator | 6. Temperature Unit |
| 3. Temperature Status | 7. Temperature Value |
| 4. Charging Indicator | 8. Record Icon |



Scan the "Register Me" QR code located on the SmartTag. If you are unable to scan the code, you may register your logger by visiting <https://register.mvebio.com> and entering the serial number found on lower left hand corner of the logger.



Download the 'Tec4App' available on both the App Store and Google Play Store.

Quick Start Guide

MVE Condition Monitoring




SmartTag continued



To start the SmartTag, press the START button and the temperature will show on the LCD. If no probe is attached, the logger will read ambient temperature.



To start recording data, press and hold the START button until the record icon  is displayed on the left side of the LCD. The SmartTag is now recording data. Log into the app or cloud to verify after registration is completed.



Stop data recording in the MVECloud or Tec4App. The STOP button is deactivated to prevent accidental stoppage during shipment.



Verify the temperature via the LCD on the SmartTag.

Once registration is completed, you will receive your login information for the MVECloud. Data is not uploaded to the MVECloud until after registration is complete.

Note: If the battery is not fully charged prior to the first shipment, please see instructions in the Operation Manual for charging the battery.

Operation Manual and additional tutorials for navigating the MVECloud and Tec4App are available at www.mvebio.com.